

Photopette[®] User Manual V1.0.2

Information in this document is subject to change without prior notice.

LEGAL DISCLAIMER

BY ACCESSING PHOTOPETTE® DEVICE AND/OR ITS APP, AND ANY SERVICES THEREOF, YOU AGREE TO BE BOUND BY THE TERMS AND CONDITIONS BELOW. THIS IS ALSO APPLICABLE FOR TERMS OF USE FOR THIRD PARTY SERVICES AND PRODUCTS INCLUDING COMPUTATIONAL PLATFORMS, REAGENTS, AND OTHER LABORATORY CONSUMABLES. BY USING ANY OF THE PRODUCT OR SERVICES OFFERED BY TIP BIOSYSTEMS PTE LTD, THE USER AGREES TO THE FOLLOWING TERMS AND CONDITIONS.

LIMITED WARRANTY

The information and materials provided in this document and related Photopette® products including text, graphics, links or other items, are provided “as is” basis. Tip Biosystems Pte Ltd is not providing any representations and warranties of any kind as to the accuracy, adequacy or completeness of such information and materials presented (or its corresponding material), and expressly disclaims liability for errors or omissions. No warranty of any kind, implied, express or statutory, including but not limited to the warranties of non-infringement of third-party rights, title, merchantability, fitness for a particular purpose and freedom from computer viruses or other harmful code, is given in conjunction with the information and materials.

COPYRIGHT

The information presented in the document and corresponding products by Tip Biosystems Pte Ltd are subject to copyright by their respective owner. Documents, content or images contained in this app, other than user generated data, may not be copied, reproduced, transmitted, displayed, distributed, altered or used in any manner without the prior written consent of Tip Biosystems Pte Ltd.

TERMS OF USE

No liability on Tip Biosystems Pte Ltd for actions related to use of product/services.

TRADEMARKS

The trademarks Photopette® and CuveTip™ are properties of Tip Biosystems Pte Ltd.

COMPANY INFORMATION

Tip Biosystems Pte Ltd (Manufacturer)

No 21, Bukit Batok Crescent, #14-79 WCEGA Tower

Singapore 658065

Tel: +65 3109 1812

Web: www.tipbiosystems.com

Declaration of conformity for the Photopette®

This is to certify that Photopette® conforms to the requirement of the following directives:

2014/35/EU	Low Voltage Equipment Safety Directive
2004/108/EC	EMI/EMC Directive
IEC 60529	Protection Class IP20

Standards to which conformity is declared, where relevant, are as follows:

IEC/EN 61010-1:2010	Safety requirements for electrical equipment for measurement, control and laboratory use. General requirements.
EN61326-1:2012	Electromagnetic compatibility- generic emission standard electrical equipment for measurement, control and laboratory use.
IEC 62133 and UN38.3	Battery certification and transport test
FCC ID: <i>PVH0950</i>	This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions; (1) This device may not cause harmful interference, and (2) This device must accept any interference received, including interference that may cause undesired operation.
IEC62471	LEDs are tested against this standard and is classified accordingly.

Subject to change and approval at the time of launch of product.

Contents

LEGAL DISCLAIMER	ii
LIMITED WARRANTY	ii
COPYRIGHT	ii
TRADEMARK	iii
Company Information	iii
Declaration of conformity for the Photopette®	iv
Introduction	7
Manual Guide	8
Photopette®	9
Device Operation	9
Accessories	11
Standard Accessories – Connecting Cable	11
Standard Accessories – Standard Solutions	11
Connectivity	11
Photopette® Technical Specification	12
Setup	15
Safety Information	15
Unpacking, Installation and Positioning	16
Safety Labels	17
Servicing Instruction	17
Application Installation	18
Sample Measurement Basics	19
Sample Handling Tips	19
Battery Operation	20
Application Usage Overview	21
Photopette® Application	21

Buttons	21
Basic Operation.....	22
Error Dialogs.....	22
Measurement Setting	23
Selecting Measurement Types	25
Selecting Datasets	25
Auto Photo Capturing.....	26
Battery Level.....	26
Using Trigger Button on Photopette Device.....	27
Measurement Screen.....	27
Datasets	29
Setting	32
Documentation and Support	33
Photopette® Applications Nucleic Acid - DNA/RNA	34
Calculations	34
Direct Protein	35
Calculations	35
Cell Viability - Resazurin/Alamar Blue™	36
Turbidity – McFarland for Cell Count.....	36
Bradford Assay	37
Other Applications.....	37
Troubleshooting.....	38
Assistance	38
Report Problem.....	38
Maintenance.....	38
Warranty	39
Cleaning and General Care.....	39

Introduction

Photopette® Photopette® is a handheld fixed-wavelength spectrophotometer for rapid, on-spot analysis. It works together with a smartphone or computing device using a dedicated application which connects via a Bluetooth™ link for taking measurements and data-transfer. It is a mobile, simple to use UV/VIS spectrophotometer instrument with a unique measurement tip (CuveTip™) coupled to a detector for accurate and precise measurement.

All Photopette® models are able to measure in small sample volume (45 µl), without any sample lost. Due to its small footprint, Photopette® can conduct tests in a very small space setting such as a fume or cell culture hood.

All calculations for a measurement are performed by a mobile device through the Photopette® App. The App is designed for using different pre-programmed Application Types with a high degree of flexibility and accessibility.

Photopette Models

509-PB – Photopette® Bio

509-PC – Photopette® Cell

509-PT – Photopette® Turbidity

509-OD600 – Photopette® OD600

509-PV – Photopette® Custom

Manual Guide

Overview This user guide describes how to operate the Photopette® device.

User documentation The guides listed below are available with the Photopette® device.

Guides
<i>Photopette® User Manual V1.0.2 (this manual)</i>
<i>Photopette® Quick Start</i>
<i>Photopette® Brochure</i>

Text and keyboard convention Text and keyboard conventions used in the *Photopette® User Manual* are listed below.


Convention	Use
Bold	Bold Text indicates user action. For example – click Run .
»	Right arrow symbol » indicates a menu choice and separates successive commands you execute or select from a drop-down or shortcut menu. For example: Select Settings » Measurement Types .

User attention words Two user attention words appear in *Photopette® User Manual* documentation. Each word implies a particular level of observation or action as described below.

NOTES: Provides information that may be of interest or help but is not critical to the use of the product.

IMPORTANT! Provides information that is necessary for proper instrument operation, accurate installation, or safe use of a chemical.

Device Operation

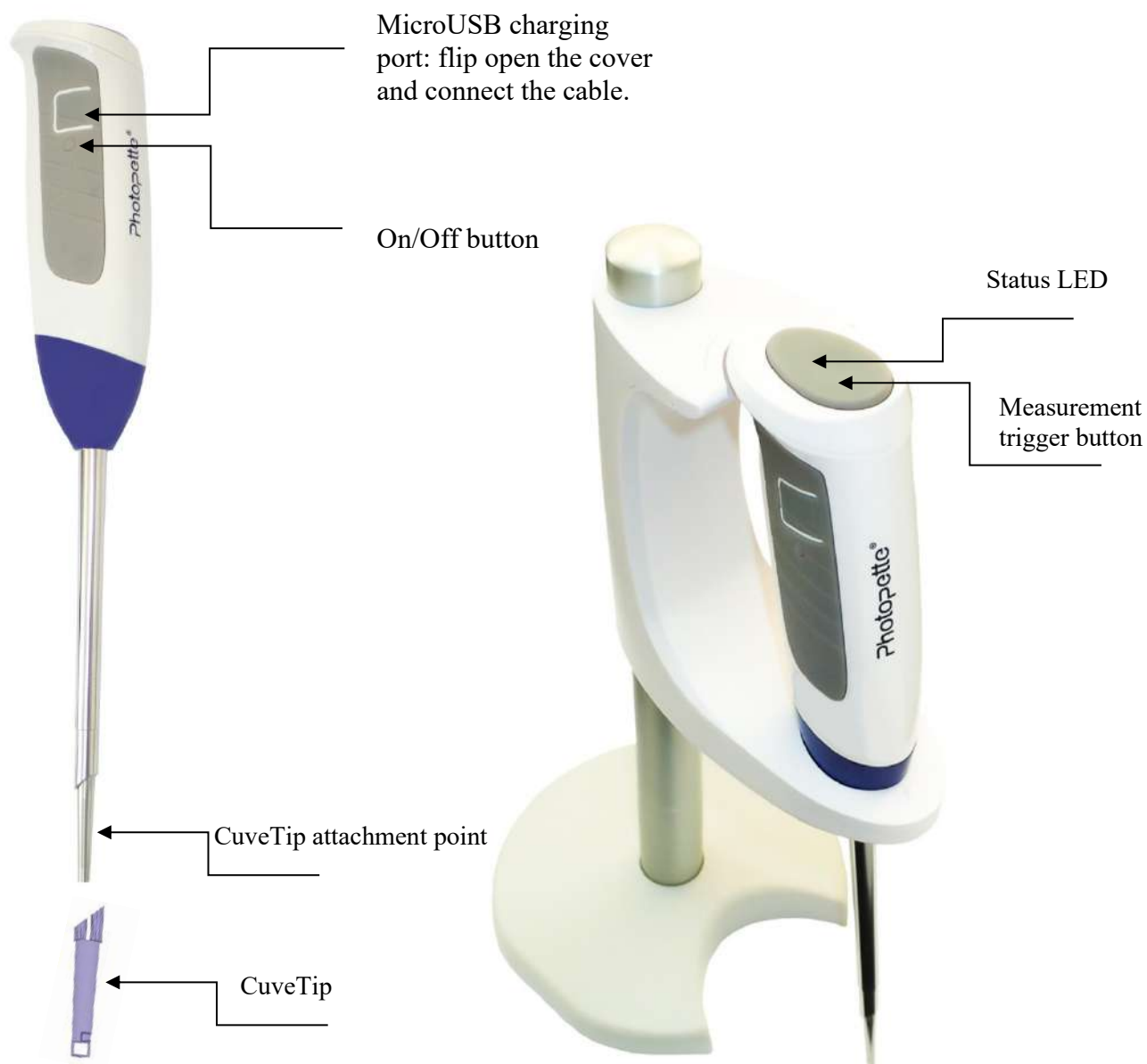

Accessories

Standard Accessories – Connecting Cable



The provided USB cable is meant for charging the Photopette® Device. Photopette® can be charged by connecting to a standard USB port capable of providing a current of 300 mA and above. A standard USB smartphone charger or a USB port on a computer/laptop is capable to supply such current.

Standard Accessories – Standard Solutions

Depending on the application type and product subscription, your Photopette® may arrive together with standard solutions to be used for quality control and maintenance purposes. **NOTES:** Proper storage and handling procedures should be followed while handling and storage of standard solution. Expired or poorly stored standard solutions should not be used for testing or quality assurance purposes.

IMPORTANT! Please read the Material Safety Data Sheet carefully prior to using this product.

Connectivity

Bluetooth

Photopette® connects to smart devices using Bluetooth™ BLE 4.0 using iOS™ and Android™ apps.

Photopette® Technical Specification

Device Specification

Interface/Connectivity	Bluetooth® BLE 4.0
Technology	Handheld fixed-wavelengths LED spectrophotometer with Ambient Light Compensation System
Measurement Time	Within 2 seconds per measured wavelength.
Warm up time	None
Lifetime (light-source)	More than 500,000 measurements per wavelength

Photopette® Bio

Wavelengths:

260 nm; Accuracy = ± 2 nm; Photometric Range = 0.00 A ~ 3.00 A
 280 nm; Accuracy = ± 2 nm; Photometric Range = 0.00 A ~ 3.00 A
 340 nm; Accuracy = ± 2 nm; Photometric Range = 0.00 A ~ 3.00 A

Photopette® Cell

Wavelengths:

340 nm; Accuracy = ± 2 nm; Photometric Range = 0.00 A ~ 3.00 A
 570 nm; Accuracy = ± 2 nm; Photometric Range = 0.00 A ~ 3.00 A
 600 nm; Accuracy = ± 2 nm; Photometric Range = 0.00 A ~ 3.00 A

Photopette® Turbidity

Wavelengths:

570 nm; Accuracy = ± 2 nm; Photometric Range = 0.00 A ~ 3.00 A
 850 nm; Accuracy = ± 2 nm; Photometric Range = 0.00 A ~ 3.00 A

Photopette® OD600

Wavelengths:

600 nm; Accuracy = ± 2 nm; Photometric Range = 0.00 A ~ 3.00 A

Photometric Precision ± 0.03 A at 1 A

Photometric Resolution 0.01 A

Power Rating	Battery Model: UR18650F-TBS Battery Type: Lithium Ion Battery Rating: DC 5 V (Charging), DC 3.7 V (Internally Powered), 2500 mAh 25,000 measurements and/or 6 days standby on a single charge Auto power-saver mode when not in use
Dimensions and Weight	Fits within 202 mm x 47 mm x 42 mm; Weight = 160 g approx. Easy Sterilization with Ethanol wipes.
Sterilization and Cleaning	Water and dust resistant housing. Resistant to mild acids and solvents.

Environmental Specification

Recommended Operating Condition	22 °C at 30% RH
Operating Temperature range	0 to 50 °C at 70% RH (non-condensing)
Operating Humidity range	30 to 80% RH (non-condensing)
Storage/Transport Temperature/ Humidity range	-20 to 50 °C at 30 to 80% RH (non-condensing)

CuveTip™

Product Variations	Sterile CuveTips™; Standard CuveTips™;
Storage Conditions	Cool, dry place with storage temperature of -20 °C to 50 °C
Analyte Temp. (max)	60 °C for 10 sec
Chemical Compatibility	Soap solution + Hydrochloric acid 36% + Sulphuric acid 40% + Acetic acid > 99% + Nitric acid 65% + Caustic soda solution 50% + Ammonia solution 33% +

Methanol	+
Ethanol	+
Isopropanol	+
Acetone	+
Butanone	+
Benzaldehyde	o
Methylene chloride	+
Plant oils	+
n-Pentane	-
Heptane	-
Toluene	-
Hexane	-
Naphtha	-
Oleic acid	-

+ Resistant; o Limited-resistance; - Non-resistant

Shelf-life

1 year for Sterile CuveTips™

2 years for Standard CuveTips™

Disposal



To be recycled as other plastics

Features and specifications are subject to change without notice.

Setup

Safety Information

Before commencing installation of the unit, please familiarize yourself with setup and operating procedures highlighted in this product guide. It is highly recommended that you go through all the warning labels and symbols on your instrument and their meaning. Please read the complete user manual prior to use.

Do not open the device as this can expose the operator to UV radiations and may damage delicate fiber optics.

IMPORTANT! Risk of UV light exposure device is used without CuveTips™.

Do not use damaged cables, accessories and another peripheral with your Photopette®.

Do not expose Photopette® to strong magnetic, electrical fields, water, chemicals or any type of liquid as heavy rain or moisture.

Do not put the device into fire, as it may swell or explode (battery). Do not store or use near any type of heat source, especially temperature above 60 °C or in an explosive atmosphere.

The device should be switched-off during transportation. The on/off button has to be protected during transportation from turning on itself caused by shock or vibration.

Do not spill any biological samples on instrument components. If spill occurs, disinfect the instrument immediately following the laboratory protocols and the cleaning instruction of the instrument.

Unpacking, Installation and Positioning

Check the contents of the package against delivery note. If any shortages are discovered, inform your supplier immediately.

Inspect the instrument for any signs of damage caused in transit. If any damage is discovered, inform your supplier immediately.

Ensure your installation site conforms to the environmental conditions for safe operation: indoor use or outdoor environment.

NOTE: Do not expose your Photopette® near liquids, chemicals, rain, moisture or dusty environments.

If the device is subjected to extreme temperature changes, it may necessary to allow the instrument to equilibrate. Turn the device off and then on again once thermal equilibrium has been established (~2-3 hours).

Avoid direct sunlight as it may bleach parts of the instrument and can cause damage plastic parts.

Please contact original supplier immediately if technical or sample handling difficulties are experienced.

NOTE: The instrument's and the accessories warranty will be void if this equipment is used in a manner not specified, or in environmental condition not suitable for safe operation. It may also adversely influence the working of the system.





Photopette Stand Installation

The Photopette stand is designed of 3 parts, the baseplate, the metal stand bar and the Photopette holder. To install, screw the metal stand bar into the base plate, then slide the Photopette holder onto the metal stand bar.

The Photopette holder is height adjustable and has a lower stop position.

The Photopette stand base plate can be connected to the CuveTip Box and the tip disposal box with the supplied clamps.

Safety Labels

Symbols	Name	Description
	UV Light caution (RG1)	Do not look directly into the distal part of the tip. This is to prevent contact with the UV light.
	Read Manual before use	User should read the user manual before use of the device so as to be aware of the precautions and proper handling of the device.
	Do not dispose in bin	Electronics are not to be disposed in the bin, this could harm the environment. Devices are to be disposed properly into electronic recycling bin at designated locations.
	Caution sign	To alert user to the relevant safety labels on the device.

Safety labels are subjected to addition and change without any notice.

Servicing Instruction

There is no servicing required for Photopette® device. Refer to user manual before use and full demo and training will be done upon delivery of devices.

Risk that might affect service personnel May get exposed to UV light from the device

Protective measures for above risk Proper service training and necessary safety labels

Verification of safe state Not required

Application Installation

The Photopette® Application is available for free download in the Play Store® and App Store. Search for “Photopette” to find the app.

Platform and compatibility	<ol style="list-style-type: none"> 1. iPhone®, iPad® or iPod® with minimum iOS 9. 2. Android phone and tablets with minimum Android 4.3 3. Photopette Application https://play.google.com/store/apps/details?id=com.tipbiosystems.noellay.photopette https://itunes.apple.com/app/id1161616983
Connectivity	<ul style="list-style-type: none"> • Internet connection with ability to download from Play Store® and App Store®.
NOTES:	<ul style="list-style-type: none"> • Photopette® Application can be downloaded through App Store®. • User will log in using the credentials to be a registered user. This allows user to be eligible for warranty coverage for device. • User will receive a confirmation email upon signing up and a notice of warranty highlighting the expiry date will be provided in the email. • Subsequently, user can start using the Application for their measurements.

Please ensure there is an established internet connection on the mobile phone before downloading the application.

The device can only run using our Application available on App Store®. It is not compatible with other applications.

Measurement data is stored locally in the smart-devices’ database and can be exported to .csv file format. It is highly recommended to frequently backup the data from the Photopette® application.

Sample Measurement Workflow

1. Switch on Photopette.
2. Start the Photopette app; enable Bluetooth.
3. Under “Device” connect with Photopette.
4. Under “Measurement type” select “Wavelength Type” or “Application Type”. Use vortexer to mix your sample to achieve homogeneous sample.
5. Attach the CuveTip onto the tip of the device. Ensure it sits tightly on the stem.
6. Press “START MEASUREMENT”.
7. Insert the CuveTip into your blank solution or blank buffer or distilled water. Make sure no bubble stays in the cavity of the CuveTip. Please read our Technical Note “How to use the CuveTip Correctly” available at the Tip Biosystems webpage at www.tipbiosystems.com
8. Press “AUTO ZERO” on the app or press the top button on the Photopette device.
9. Remove CuveTip from the blank. In case you sample has a large volume of more than 10 mL you can directly insert the same CuveTip into your sample and move the tip for 10 seconds to exchange any blank solution residues with sample. For small sample volumes, tap the CuveTip onto a clean lab paper to remove the blank by capillary force first before transferring into the sample.
10. Press “MEASUREMENT” and the absorbance values or concentrations will be shown directly on the screen.
11. You can use the same CuveTip to perform additional measurements.

NOTE: Do not overfill the sample vial. This is to prevent samples from seeping in from the top of the sample vial.

NOTE: Check the chemical compatibility chart for the compatibility of the sample solvent or solution you are using. The information is available in the annexure of this manual and on the technical information of the website. Please contact us at support@tipbiosystems.com if your solvent of your sample is not listed on the chart.

Sample Handling Tips

- It is important to mix or vortex every sample right before measurement, if needed, to ensure homogeneity.
- The minimum volume that can be used for sample measurement is 45 µl and can be reuse if it is inserted into the tip cavity using a micropipette.
- Proper fitting of the CuveTip™ is important to ensure accurate measurement.
- In most cases, a laboratory clean wipe is sufficient to clean the device.
- Please avoid touching the tip of the device.

Battery Operation

The device runs on rechargeable lithium-ion batteries. The batteries can be operated for approximately 12 hours within a normal usage i.e. 20-30 measurement per hour. The device has a charging indicator LED at the top button. The charging indicator will change to green once it is fully charged. When the battery gets low there will be an alert on the application to charge. Once it is below 10% there will be a beeping sound. At this status, it is still possible to use the device for another one hour. However, it is recommended to charge the battery at this status. If the battery is not charged there will be a second warning message when the battery is empty and several beeping sounds before the instrument is automatically shut down within 10 seconds.

The battery is standard Lithium-ion battery with a lifespan of about 500 full charging cycles. Subsequently, the capacity and operation time of the battery may vary. A fully charge battery will self-discharge over a time of 30-40 days. After this time recharging of the battery is necessary. Charging time for an empty battery is 5 hours.

NOTE: If Photopette shuts down during a measurement process, the data of the last measurement will be still retained in the dataset.

Application Usage Overview

Photopette® Application

Main Features	Measurement
	Setting
	Measurement
	Dataset page
	Setting

NOTE: User should set **Measurement Setting** before starting measurements. The user can view their measurements by dataset names on the **Datasets page**. In the Datasets page, user can also rename datasets, delete datasets, and share datasets. User can test device performance, change languages (currently English, Chinese and German only) and read “About” in the **Setting page**.

The Photopette® App has a set of pre-programmed applications according to the type of wavelengths present in the device. To select an Application type, go to measurement type to see a list of Application types user can use.

Buttons



Prior to any sample measurement at the start of a new method, an auto zero measurement of either water or the buffer of the samples is required to give the Photopette a reference of what zero should be.

To initiate the scan of the sample, press the measurement button, the data will be stored in the Dataset until the method is exited at this time the user is able to share or delete the existing datasets.

Basic Operation

Error Dialogs

Error #	What does it mean?	What can I do?
A1, M1	The detected signal exceeds the saturation limit of the detector. This error may appear when measuring in very bright environments.	Move to a location with reduced ambient light or try to block ambient light before retaking the measurement.
A2, M2	Photopette seems to experience some technical difficulties.	Restart the device and perform the Self-Test. If the problem is still present, please contact your local distributor for support.
A3, M3	Photopette did not detect any CuveTip placed on the device. This error may also occur if the measurement is performed on a highly absorbing sample.	Ensure the CuveTip is placed properly on the device. When measuring on a highly absorbing sample, this warning can be ignored.
A4	Photopette automatically compensates for ambient light up to a certain threshold. Close to the threshold, accuracy of the auto-zero might be reduced as fewer data points are available for the compensation. You can still choose to continue your measurements.	Move to a location with reduced ambient light or try to block ambient light before retaking the measurement.
A5	Photopette automatically compensates for ambient light up to a certain threshold. It seems there are fewer data points available for the compensation than usual. You can still choose to continue your measurements.	Please verify device performance via the Self-Test. If the error re-occurs, try changing to a location with reduced ambient light.
A6, M4	Photopette did not detect any CuveTip placed on the device. This error may also occur if the measurement is performed on a highly absorbing sample. You can still choose to continue your measurements.	Ensure the CuveTip is placed properly on the device. When measuring on a highly absorbing sample, this warning can be ignored.


Measurement Setting

User should set Measurement Setting before measurement process. This allows the device to store the required settings for the measurements.

Measurement Settings:

- Selecting Device
- Selecting Measurement Type
- Selecting Dataset
- Auto Photo Capturing
- Auto Measurement
- Battery Level

Selecting Device

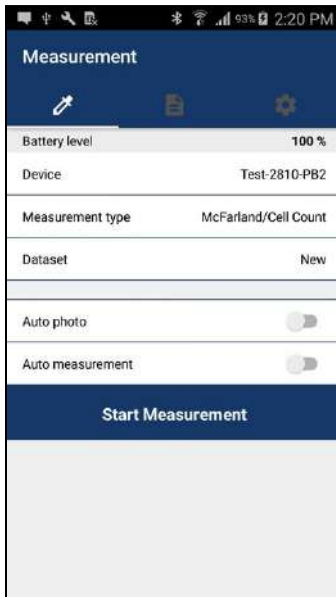

Before start measurement, most important thing to do is to connect with the Photopette® device.



Users must **ON** Bluetooth connection in their device in order to connect with the Photopette® device.

1. Connecting with Photopette™ Device (Figure 4.1) from the Application

- a. **Measurement Setting » Device**
- b. Available Photopette® devices is listed
Select the correct Device listed.
- c. If connection successful, device will make 3 beeping sound and note the tick mark on right side of connected device name.

2. Renaming Photopette® Device

- a. User can rename photopette device by swiping device name view.
- b. By right swiping device name view, it will come out **Edit** button and **Disconnect** button. (Disconnect button is only available for connected device.)
- c. Please click **Edit** button and fill in proper name at “Rename” window.
- d. Click **OK** and user can see device name already change.
- e. For subsequent connection, the user can see a personalised name on their device. This will help in identifying and differentiating multiple devices.

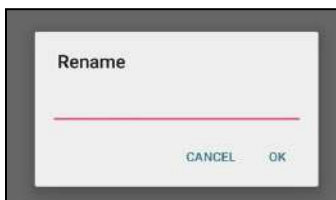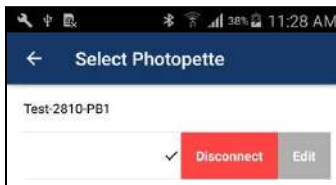

Selecting Measurement Types



In Measurement Setting » Measurement Type

Select Option -Wavelength Type

-Application Type

NOTES:


If user choose one application type, one or more wavelength that corresponding to this application type will be measured.

For instance,

If user chooses ssDNA, 260 nm and 280 nm wavelength will be measured.

According to Measurement type that user choose, selectable datasets will be different.

Selecting Datasets

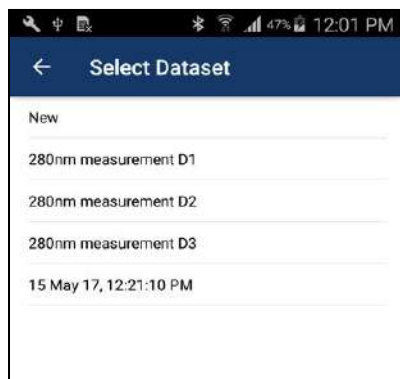

To store measurement information, the user needs to select a dataset first.

Select Measurement Setting » Dataset

Select Option - New
 - Existing Datasets

NOTES:

If user select **New Dataset**, new dataset will be created to save measurement information.

If user select one of existing dataset, information will be appending on information of selected dataset.

User can see existing datasets with same measurement type that user chooses.

Auto Photo Capturing



If user would like to automatically capture photo on measurement time, **Auto photo** should be enabled.

Auto Measurement



If user would like to automatically measure, please enable **Measurement Setting » Auto measurement** (Figure 4.3)

User can see 3 additional settings by enabling Auto measurement.

1. Numbers of Reading
2. Time Interval
3. Alert me when done

NOTES: Please choose appropriate time interval depending on number of wavelengths to be measured and Auto photo.

Time Interval = No of wavelength *2 + Auto Photo

For instance, if user chooses dsDNA, 260 nm and 280 nm will be measured. So, time interval is 4sec. If auto photo is On, application will need 3 more sec to take photo automatically. So, final time interval needed is 7 sec.

Please check battery level often. If battery is lower than 40%, please recharge in order to ensure proper measurement.

Battery Level

Using Trigger Button on Photopette Device

Activating Trigger Button

Instructions

1. Click once in Setting to start measurement
2. Click for auto zero and you can hear beep sound of measurement event.
3. Click for warning/dialog dismissal.
4. Click for final measurement.
5. User can view Transmittance and Absorbance value on measurement screen.
6. For additional measurement, click once again.

IMPORTANT: Please click and listen beep sound of recognition. Continuous clicking will result in many measurements.


Measurement Screen

Measurement Screen Functions

1. Taking Photo
2. Adding Note
3. New Measurement
4. Cancel Auto Measurement
5. View Measurement Results

Components

1. Measurement Count
2. Last measurement time
3. Photo/Notes Button
4. Measurement/ Cancel Button
5. Measurement Result Table
6. Back Button
7. Dataset button

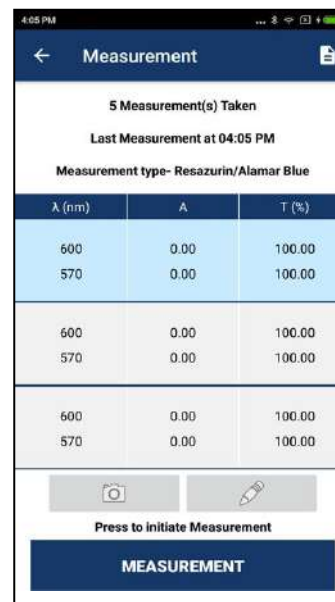

1. Taking Photo Click **Photo**  and capture photo before measurement start.

NOTE: If user enable **Auto Photo** or **Auto Measurement** in Setting, **Photo** button in measurement screen will be disabled.

2. Adding Note Click **Notes**  and add note before measurement start.

NOTE: If user enables **Auto Measurement** in Settings, **Notes** button in measurement screen will be disabled.

3. New Measurement User can repeat measurement manually by clicking Measurement after a measurement finished.

NOTE: If user enable **Auto Measurement** in Setting, **Measurement** button in measurement screen will not be seen.


4. Cancel Auto Measurement User can cancel auto measurement anytime by clicking **Cancel** on Measurement Screen.

5. View Measurement Results Transmittance and Absorbance result will be shown in every measurement.

Concentration result will be shown for some application type only.

Warning icon will be shown for measurement accuracy.

User can click Warning icon and see what is warning.

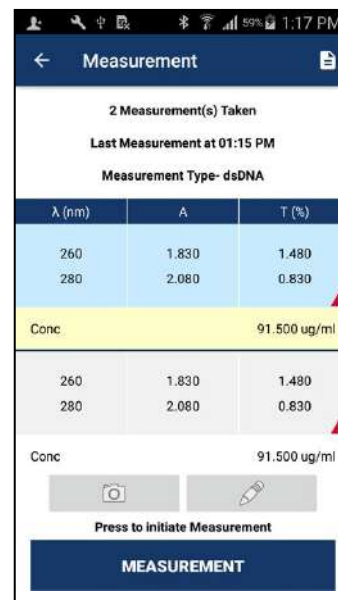

Datasets

All measurement result will be stored in a dataset.

Dataset Functions

1. View Dataset List
2. View Detail Measurement
3. Rename Dataset
4. Share/Delete Datasets

1. View Dataset List

Select **Datasets** from navigation bar.
Datasets list that user measured will be shown.


NOTE: Auto Dataset name is Measurement Time DD MMM YY, hh:mm:ss AM/PM.

2. View Measurement Details

Select **Datasets** details by clicking


Measurement Detail Information:

- Measurement Date and Time
- Measurement Type
- Wavelength
- Absorbance value
- Transmittance
- Concentration (if available)
- Photo (Enable if photo is available)
- Notes (Enable if note is available)
- Warning (if available)



Measurement type - Resazurin/Alamar Blue			
λ (nm)	A	T (%)	
27 Feb 17, 04:05:12 PM			
600	0.00	100.00	
570	0.00	100.00	
27 Feb 17, 04:05:06 PM			
600	0.00	100.00	
570	0.00	100.00	
27 Feb 17, 04:05:00 PM			
600	0.00	100.00	
570	0.00	100.00	
27 Feb 17, 04:04:44 PM			
600	0.00	100.00	


3. Rename Dataset

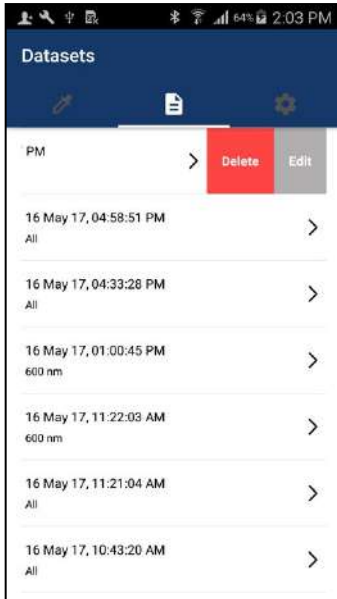

1. User can rename dataset by swiping a dataset view.
2. By right swiping a dataset view, **Delete** button and **Edit** button will appear.
3. Please click **Edit** button and fill in proper name at “Rename” window.
4. Click **OK** and user can see the change of the device name.

After rename is completed, the user can see a customised dataset name. This will help in identifying and differentiating multiple datasets.

4. Share/ Delete datasets

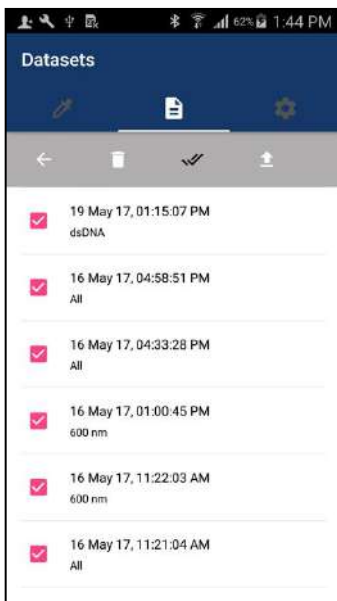

User can share/delete datasets with 2 options:

1. Share/Delete All Datasets
2. Share/Delete two or more datasets

Android



1. Share/Delete All Datasets

Long press a dataset to appear toolbar

Click **Tool bar** » **Select All** 

For Share datasets, Click **Tool bar** » **Share** 

Two Share Options will appear:

1. Email 
2. Google Drive 

For Delete datasets, Click **Tool bar** » **Delete** 


2. Share/Delete one or more datasets

Long press a dataset to appear checkbox

Select dataset item that should be deleted or shared.

Click **Tool bar** » **Share or Delete**


Selected datasets can be shared or delete.

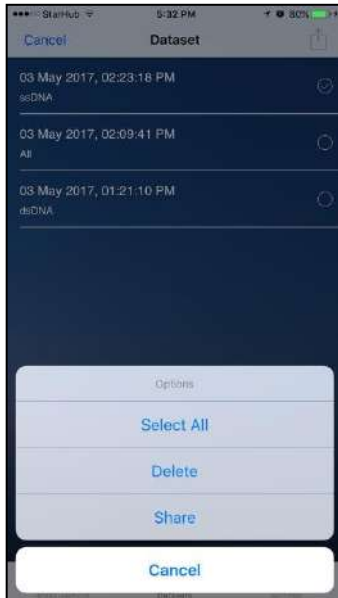


iOS

3. Share/Delete All Datasets


Long press a dataset or click  to appear check boxes

Click , user will see Select all, Delete and Share options.

For Share datasets, Click  » **Share**

Two Share Options will appear:

- 3. Email
- 4. Google Drive

For Delete datasets, Click  » **Delete**

4. Share/Delete one or more datasets

Long press a dataset to appear checkbox

Select dataset item that should be deleted or shared.

Click  » **Share or Delete**

Selected datasets can be shared or delete.

NOTE: Dataset CSV format file will be exported to google drive/email.

If user would like to deselect all from **Select All**, just click **Deselect** on top bar.

Setting


Setting page offers additional features that does not connected with measurement process.

Setting

User can perform 3 features in Setting.

1. Self-Test
2. Change Language
3. About

Self-Test

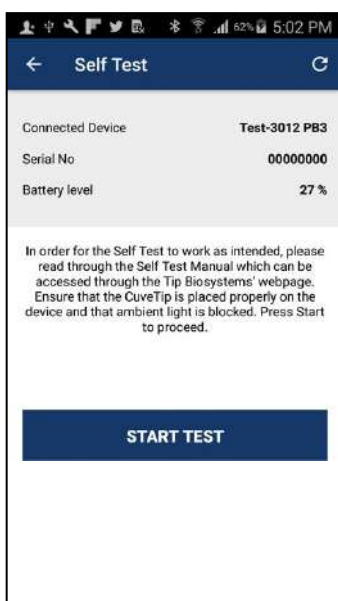

Self-Test allow users to test device performance by themselves. User can follow step-by-step guideline that prompt on app and can get device condition as a final result.

The following information are shown default in Self-Test:

1. Connected Device Name
2. Serial Number
3. Battery Level

NOTE: Self-Test can only perform when app is connected

Change Language



User can change 3 languages in Change Language.

1. English (Default)
2. Simplified Chinese
3. German

NOTE: By changing language, all text within the app will show according to changed language.

About



1. App Version
2. Firmware Version
3. Date of Release
4. Legal Disclaimer
5. Limited Warranty
6. Copyright
7. Limitation of Liability

NOTE: Item 2. Firmware Version is only shown when device is connected with Photopette app.

To view Item 4,5,6,7 Select About » Read Legal Disclaimer

Documentation and Support

Obtaining Support

Technical support

For any technical questions, please contact us via email – support@tipbiosystems.com

Limited Product Warranty

Tip Biosystems Pte Ltd and/or its affiliate(s) warrant their products as set forth in the Tip Biosystems' General Terms and Conditions found on Photopette Application About Page.

If you want to know more, please contact Tip Biosystems Pte Ltd at <http://www.tipbiosystems.com>

Photopette® Application Examples

A full list of all applications can be found on the Tip Biosystems webpage at: www.tipbiosystems.com

Nucleic Acid - DNA/RNA

Measurement of DNA and RNA concentration and determination of DNA and RNA purity can be done by the measurement of absorbance at the wavelength of 260 nm and 280 nm.

In the iOS application, under the **Measurement Type**, select the nucleic acid category (DNA or RNA).

Calculations

Example:

A DNA sample has an absorbance reading of 0.4. The concentration is calculated as:
 $0.4 * 50 = 2 \mu\text{g/ml}$

Nucleic Acids	Concentration ($\mu\text{g/ml}$) per A_{260} Unit
dsDNA	50
ssDNA	33
ssRNA	40

Table 1: Calculations for nucleic acid concentration

Purity of	A_{260}/A_{280} Ratio
DNA	~ 1.8 (approximate value)
RNA	~ 2.0 (approximate value)

Table 2: Calculations for determination of nucleic acid purity

For purity check of nucleic acid, a ratio of A_{260}/A_{280} is used.

A ratio lower than 1.8 for DNA and a ratio lower than 2.0 for RNA indicates presence of contamination. However, the A_{260}/A_{280} ratio is only an indication of purity rather than a precise answer.

NOTE: Small changes of the pH of the solution will affect the 260/280 ratio. Acidic solution will under represent the 260/260 ratio (0.2-0.3) and basic solution will over represent the ratio.

NOTE: It is important to check for the pH of a sample measured on our instrument as acidic and basic solutions will vary the absorbance value.

Direct Protein

In the iOS application, under the Measurement Type, select the Direct Protein.

Total protein concentration can be determined by the measurement of absorbance at 280 nm. This is a simple and quick way of quantification and the protein sample is recoverable. Proteins absorb light at 280 nm, primarily the aromatic amino acids (tryptophan and tyrosine), and the disulfide bonds (S-S bonds) in the protein sequence. The amount of light absorbed by the protein is dependable on the chemical composition of the protein – the number and type of amino acids. Protein absorption is also affected by the protein structure. Therefore, conditions such as temperature, pH ionic strength, and presence of detergents can affect the absorption of the proteins.

Calculations

To compute the concentration of a purified protein solution, the formulas are given as:

1. A_{260}/A_{280} Ratio: To check for a pure protein, the A_{260}/A_{280} Ratio is approximately about 0.58.
2. To eliminate the presence of nucleic acids in the protein sample, the protein concentration can be estimated according to Warburg and Christian formula:

$$\text{Concentration (mg/ml)} = (1.55 \times A_{280}) - (0.76 \times A_{260})$$

3. For better calculation of proteins:

$$\text{Protein concentration (mg protein or peptide per ml)} = (A_{280} \times \text{DF} \times \text{MW}) / \epsilon$$

Where A_{280} is the absorbance of the solution at 280 nm in a 1-cm cell

DF is the dilution factor

MW is the molecular weight of the protein/peptide

ϵ is the molar extinction coefficient of each chromophore at 280 nm

Cell Viability - Resazurin/Alamar Blue™

Tip Biosystems offers the measurement of the reduction of resazurin to resofurin. In this cell viability assay, the resazurin exhibits an absorption peak at 600 nm and the resofurin at 570 nm wavelengths. An alternate wavelength measurement for absorption peak is available at 630 nm.

Resazurin is an indicator dye which incorporates an oxidation-reduction (REDOX) indicator that changes color in response to the chemical reduction of growth medium of the growing cells. The bioassay can be used to establish proliferation or relative cytotoxicity. The toxicity of related novel agents can be predicted by comparing its baseline data against the baseline data of the known *in-vivo* toxicity.

Turbidity is the cloudiness or haziness of a fluid caused by large numbers of individual particles that are present but undissolved. Determining Turbidity of an aqueous solution is a frequently used procedure to estimate the quality of water. Turbidity is determined by measuring the optical density (OD) at 600 nm since light is scattered rather than absorbed. Careful mixing of the suspension is required before each measurement is taken.

Turbidity - A calibration graph can be plotted using Turbidity 4000 NTU Calibration Standard – Formazin from Sigma Aldrich (Product number TURB4000). The turbidity of the test solution is then determined by substituting the OD value in the equation of the straight line that is determined from the calibration graph.

Turbidity – McFarland for Cell Count

The number of cells in a sample can be estimated by measuring the OD of the cells suspended in a liquid. A calibration curve from the absorbance of known cell density is plotted and the absorbance of the unknown cell density is compared against the calibration curve. This is a quick way to determine the number of bacteria in a sample. The OD value may differ when different type of cells is used. Therefore, the calibration curve can only be used for the same type of cell.

McFarland Standards are usually labeled as 0.5 to 10. They can be made in the lab by preparing a 1% anhydrous BaCl₂ and 1% solution of H₂SO₄ and mix both solutions at specified volumes and ratios. Commercially available latex beads suspensions are also available.

A calibration graph is plotted from the OD of the McFarland Standards. The turbidity of the test solution is then determined by substituting the OD value in the equation of the straight line that is determined from the calibration graph.

Bradford Assay

In the iOS application, under the Measurement Type, select the Bradford Assay. The Bradford Assay is used to determine protein concentrations. The Bradford reagent contains the Brilliant Blue G dye which will form a colored complex with the protein in the solution. The linear concentration range is 0.1 – 1.4 mg/ml of protein, using BSA (bovine serum albumin) as the standard protein and Bradford reagent (Product number B6916) from Sigma-Aldrich. A standard curve is generated using serial dilutions of known BSA concentrations. An unknown concentration of a protein can be determined by comparing against the standard curve.

Other Applications

Other applications are available as Application Notes for downloading at Tip Biosystems Pte Ltd webpage www.tipbiosystems.com. For applications not listed in this manual or at Tip Biosystems's webpage please contact our application specialist at support@tipbiosystems.com

Troubleshooting

The Photopette App has a self-test function under “Settings”. The Photopette® calibration test is performed with the standard solutions provided in the package. If the instrument passes the self-calibration test, the application will reflect the outcome. If the device does not pass the test, an alert will appear on the application along with the message explaining the reason for the failed test along with the recommended solution.

If the Ok button is selected, the window is closed and the home screen will be shown with an error message at the top. The error message is also shown on top of all other screen and all performed measurements including the saved files of those measurements. If the self-calibration test fails, please contact the support@tipbiosystems.com.

Assistance

The assistance includes off-site support, report a problem, software maintenance and legal as function to help with any technical issues or questions that may arise with the Photopette®.

Report Problem

The Application will send bug error reporting automatically to our Software Developer in case there is any problem or the application crashes. Other related problems can be emailed to the above email accordingly.

Maintenance

Download the software update file from Tip Biosystems main download page. Remember to save all data before updating the application. Please ensure that your mobile device has enough storage and is connected to Wi-Fi before commencing Application download.

Warranty

Tip Biosystems Pte Ltd guarantees that the product supplied has been thoroughly tested to ensure that it meets its published specification. The warranty included in the conditions of supply is valid for 12 months and is valid only if the product has been used according to the instructions supplied.

Tip Biosystems Pte Ltd or your supplier can accept no liability of loss or damage arising from the faulty or incorrect use of this product. Each device will come with a Certificate of Performance where upon assembly.

Cleaning and General Care

Switch off the Photopette® and disconnect the power cord prior to external cleaning. Use a soft wet cloth or dry microfiber cloth to clean all external surfaces. A mild liquid detergent may be used to remove stubborn marks.

The Photopette device can be sterilized by ethanol solution applied with a wet cloth. Do not over apply.

Approved disinfectant solutions include: Apesin disinfection spray (Tana Chemi GmbH), Incidin Liquid & Inciddin Foam (Ecolab), and Lysoformin Spezial (Lysoform Dr. Hans Roseman GmbH).

NOTE: Observe all necessary precautions if dealing with hazardous samples or solvents.